

PCR protokoll

Datum: 120316

Projekt: Acoela 2011

Primers: LCO/HCO



Program:

	ul/prov	mix X
H ₂ O	22	
primer 1	0,5	
primer 2	0,5	
DNA	2	

95°C	4min	X1
94°C	30"	X44
46 °C	30"	
72°C	30"	
72°C	10 min	X1

Maskin nr:

1	UJ11_20_	26	UJ11_314_	51	UJ11_393_
2	UJ11_21_	27	UJ11_315_	52	UJ11_398_
3	UJ11_35_	28	UJ11_317_	53	UJ11_225II_
4	UJ11_36_	29	UJ11_318_	54	kontroll
5	UJ11_72_	30	UJ11_319_		
6	UJ11_75_	31	UJ11_320_		
7	UJ11_76_	32	UJ11_324_		
8	UJ11_79_	33	UJ11_325_		
9	UJ11_82_	34	UJ11_329_ <i>gds</i>		
10	UJ11_111_	35	UJ11_330_		
11	UJ11_155_	36	UJ11_331_		
12	UJ11_159_	37	UJ11_332_		
13	UJ11_167_	38	UJ11_333_		
14	UJ11_172_	39	UJ11_341_		
15	UJ11_202_	40	UJ11_344_		
16	UJ11_211_ <i>gds</i>	41	UJ11_356_		
17	UJ11_213_ <i>gds</i>	42	UJ11_357_		
18	UJ11_267_	43	UJ11_362_		
19	UJ11_272_	44	UJ11_363_		
20	UJ11_276_	45	UJ11_365_		
21	UJ11_277_	46	UJ11_379_ <i>gds</i>		
22	UJ11_302_	47	UJ11_382_		
23	UJ11_303_	48	UJ11_386_		
24	UJ11_305_	49	UJ11_387_		
25	UJ11_309_	50	UJ11_388_		

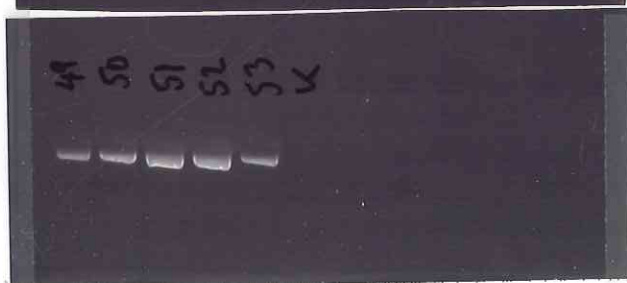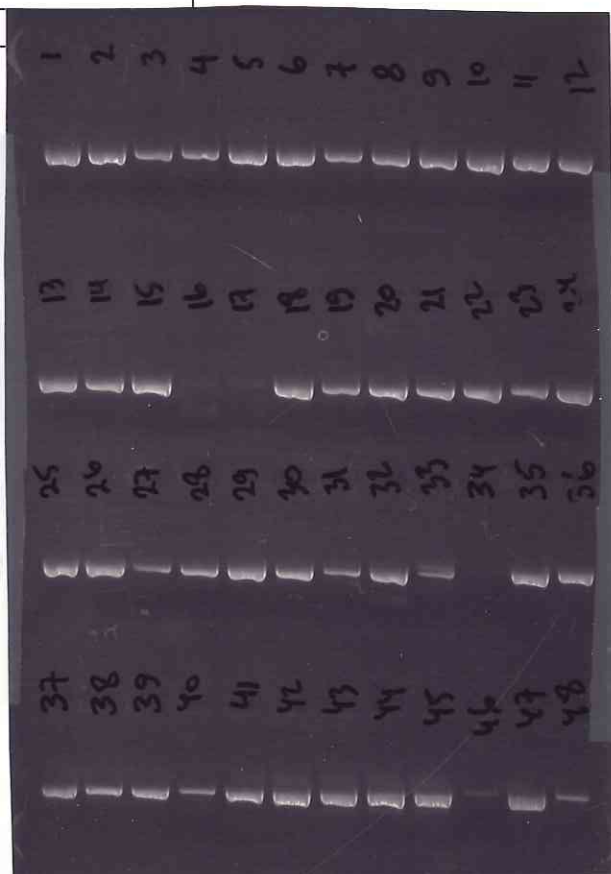

Keyvan

* Om PCR på alla extrakt
med nybeställda primer; 120112
Kontroll neg